



PRODUCT DESCRIPTION

AS-1267 | LB Agar (Miller)
www.ausamics.com.au

LB Agar (Miller)

High-Salt Bacterial Culture Medium for Routine Molecular Biology

Catalogue Number: AS-1267

Application	Medium Type	Typical pH	HS Code
E. coli cloning & routine culture	Non-selective solid agar medium	7.0 ± 0.2 at 25 °C	3821.00.00

Overview

LB Agar (Miller) is the most widely used general-purpose solid culture medium for the routine cultivation, cloning, and maintenance of recombinant *Escherichia coli* strains in molecular biology laboratories. The Miller formulation, originally described by J.H. Miller in 1972, contains 10.0 g/L sodium chloride — higher than the Lennox (5.0 g/L) and Bertani (0 g/L) variants — making it the global standard for routine *E. coli* work.

This formulation is fully compatible with common antibiotics, supplements, IPTG/X-gal reporter systems, and a wide range of laboratory strains. It supports rapid bacterial growth, high cell density culture, and reliable plasmid maintenance.

Principle of the Medium

Tryptone (10.0 g/L)	Pancreatic digest of casein providing peptides, amino acids, nitrogen, and carbon for rapid <i>E. coli</i> growth
Yeast Extract (5.0 g/L)	Supplies B-vitamins, trace elements, nucleotides, coenzymes, and growth factors essential for <i>E. coli</i> K-12
Sodium Chloride (10.0 g/L)	High-salt Miller formulation — maintains optimal osmotic strength for standard <i>E. coli</i> strains
Agar (15.0 g/L)	Solidifying agent providing a firm, clear growth surface suitable for colony isolation and selection

Typical Composition (per litre)

Component	CAS Number	Function	Amount
Tryptone (Pancreatic Digest of Casein)	73049-73-7	Nitrogen, carbon, amino acids	10.0 g



PRODUCT DESCRIPTION

AS-1267 | LB Agar (Miller)
www.ausamics.com.au

Yeast Extract	8013-01-2	Vitamins, growth factors, nucleotides	5.0 g
Sodium Chloride	7647-14-5	Osmotic balance (high-salt Miller)	10.0 g
Agar	9002-18-0	Solidifying agent	15.0 g

Total per litre: 40.0 g | Final pH: 7.0 ± 0.2 at 25 °C

Formulation Comparison — LB Variants

Component	Miller (AS-1267)	Lennox (AS-1269)	Bertani
Tryptone	10.0 g/L	10.0 g/L	10.0 g/L
Yeast Extract	5.0 g/L	5.0 g/L	5.0 g/L
NaCl	10.0 g/L	5.0 g/L	0 g/L
Agar	15.0 g/L	15.0 g/L	15.0 g/L
Total solids	40.0 g/L	35.0 g/L	30.0 g/L
Best for	Standard E. coli — global default	Salt-sensitive strains	Very low osmolarity work

Applications

Molecular Biology & Genetic Engineering

- Routine plating of E. coli strains — DH5α, BL21, TOP10, JM109, XL1-Blue
- Plasmid propagation, transformation, and colony screening
- Antibiotic selection plates (ampicillin, kanamycin, chloramphenicol, tetracycline)
- Blue–white screening when supplemented with X-gal (40 µg/mL) and IPTG (0.1 mM)
- Recombinant protein expression colony selection

Phage & Microbiology Research

- Propagation of plasmid-containing E. coli strains
- General bacteriophage and host strain studies
- Overlay plaque assays when combined with top agar



PRODUCT DESCRIPTION

AS-1267 | LB Agar (Miller)
www.ausamics.com.au

Teaching & General Laboratory Use

- Standard medium for molecular biology instruction
- Routine microbiology cultivation and streak plate techniques
- Quality control of *E. coli* competent cell preparations

Preparation Instructions

1. Suspend 40.0 g of dehydrated medium in 1 litre of demineralised or purified water.
2. Heat with frequent agitation until fully dissolved.
3. Verify and adjust pH to 7.0 ± 0.2 if required using 1 M NaOH or 1 M HCl.
4. Sterilise by autoclaving at 121 °C for 15 minutes.
5. Cool to 45–50 °C.
6. Add filter-sterilised antibiotics or supplements aseptically at this stage if required.
7. Dispense 20–25 mL per 90 mm sterile Petri dish.
8. Allow to solidify at room temperature on a level surface before use or storage.

Note: Do not overheat or re-autoclave prepared agar. Add heat-sensitive supplements (X-gal, IPTG, antibiotics) only after cooling to 45–50 °C.

Common Antibiotic Selection Concentrations

Antibiotic	Working Concentration	Stock Concentration	Storage
Ampicillin	100 µg/mL	100 mg/mL in water	-20 °C
Kanamycin	50 µg/mL	50 mg/mL in water	-20 °C
Chloramphenicol	25 µg/mL	25 mg/mL in ethanol	-20 °C
Tetracycline	12.5 µg/mL	12.5 mg/mL in ethanol	-20 °C, protect from light
Gentamicin	10 µg/mL	10 mg/mL in water	-20 °C
Spectinomycin	50 µg/mL	50 mg/mL in water	-20 °C

Note: Filter-sterilise all antibiotic stocks through a 0.22 µm membrane. Add aseptically after cooling to 45–50 °C.

Blue–White Screening Supplements



PRODUCT DESCRIPTION

AS-1267 | LB Agar (Miller)
www.ausamics.com.au

Supplement	Final Concentration	Preparation	Purpose
X-gal	40 µg/mL	20 mg/mL in DMF or DMSO	Chromogenic substrate for β-galactosidase
IPTG	0.1 mM	100 mM in sterile water	Inducer of lac operon

Spread 40 µL of X-gal stock and 4 µL of IPTG stock onto the agar surface 30 minutes before plating. Blue colonies contain lacZ insert; white colonies are recombinants.

Incubation & Typical Growth

Inoculate by surface spread or pour plate. Incubate aerobically at 37 °C for 12–18 hours. Colonies are visible within 12–16 hours.

Organism	Expected Result	Incubation
E. coli K-12 DH5α	Good growth — white/cream 1–3 mm colonies	37 °C, 12–16 h
E. coli BL21(DE3)	Good growth — suitable for expression screens	37 °C, 12–16 h
E. coli TOP10	Good growth — high-efficiency cloning host	37 °C, 12–16 h
E. coli JM109	Good growth — compatible with M13 phage work	37 °C, 12–16 h
Salmonella typhimurium ATCC 14028	Good growth	37 °C, 18–24 h

Storage & Stability

Dehydrated powder	Store at 15–30 °C in a cool, dry, tightly closed container
Protect from	Moisture, direct light, and extreme temperatures
Prepared plates (no antibiotics)	2–8 °C, inverted, up to 4 weeks
Prepared plates (with antibiotics)	2–8 °C, use within 1–2 weeks
Shelf life (dehydrated)	As per labelled expiry date
Do not use if	Contamination, colour change, or cracking of agar surface is observed



PRODUCT DESCRIPTION

AS-1267 | LB Agar (Miller)
www.ausamics.com.au

Quality & Compliance

Manufactured under controlled conditions for batch-to-batch consistency. Performance tested against ATCC reference organisms prior to release. Compatible with protocols described in Miller (1972) Experiments in Molecular Genetics, Sambrook & Russell (Molecular Cloning), and Current Protocols in Molecular Biology.

Reference

Miller, J.H. (1972). *Experiments in Molecular Genetics*. Cold Spring Harbor Laboratory, New York.

Customs & Trade Information

HS / AHECC Code	3821.00.00
Description	Prepared culture media for the development or maintenance of microorganisms
Country of Origin	Australia

Disclaimer

This product is manufactured and supplied by AuSaMicS Pty Ltd for laboratory and research use only. It is not intended for human or veterinary consumption, therapeutic use, or in vitro diagnostic procedures without appropriate validation. AuSaMicS Pty Ltd makes no representations or warranties, express or implied, regarding the fitness of this product for any particular purpose beyond its stated intended use. Users are solely responsible for compliance with all applicable laws, regulations, and safety requirements. AuSaMicS Pty Ltd shall not be liable for any direct, indirect, incidental, or consequential damages arising from the use or misuse of this product. All information provided is believed to be accurate at the time of publication and is subject to change without notice.