



PRODUCT DESCRIPTION

AS-1270 | LB Broth (Miller)
www.ausamics.com.au

LB Broth (Miller)

Luria-Bertani Broth, Miller Modification — General-Purpose Liquid Culture Medium

Catalogue Number: AS-1270

Application	Medium Type	Typical pH	HS Code
E. coli & molecular biology	Non-selective liquid broth	7.0 ± 0.2 at 25 °C	3821.00.00

Overview

LB Broth (Miller), Luria-Bertani Broth Miller modification, is a nutrient-rich, general-purpose liquid culture medium widely used for the cultivation and maintenance of *Escherichia coli* in molecular biology and recombinant DNA applications. It is a modification of the original Luria-Bertani formulation with a higher sodium chloride concentration (10.0 g/L), which supports osmotic balance and membrane transport across a broad range of laboratory *E. coli* strains.

The combination of tryptone and yeast extract supplies essential amino acids, peptides, vitamins, and growth factors, supporting robust growth of *E. coli* strains including K-12 derivatives with auxotrophic requirements. LB Broth (Miller) is commonly used as both a liquid culture medium and as a base for LB Agar (Miller) plate preparation.

Principle of the Medium

Tryptone (10.0 g/L)	Pancreatic digest of casein — primary source of peptides, amino acids, nitrogen, and carbon for rapid <i>E. coli</i> growth
Yeast Extract (5.0 g/L)	Supplies B-vitamins, trace elements, nucleotides, coenzymes, and growth factors essential for <i>E. coli</i> K-12
Sodium Chloride (10.0 g/L)	High-salt Miller formulation — maintains optimal osmotic strength for standard laboratory <i>E. coli</i> strains

Typical Composition (per litre)

Component	CAS Number	Function	Amount
Tryptone (Pancreatic Digest of Casein)	73049-73-7	Nitrogen, carbon, amino acids, peptides	10.0 g



PRODUCT DESCRIPTION

AS-1270 | LB Broth (Miller)
www.ausamics.com.au

Yeast Extract	8013-01-2	B-vitamins, growth factors, nucleotides	5.0 g
Sodium Chloride	7647-14-5	Osmotic balance — Miller high-salt formulation	10.0 g

Total per litre: 25.0 g | Final pH: 7.0 ± 0.2 at 25 °C

Comparison: LB Broth Formulations

Component	Miller (AS-1270)	Lennox (AS-1269)	Bertani
Tryptone	10.0 g/L	10.0 g/L	10.0 g/L
Yeast Extract	5.0 g/L	5.0 g/L	5.0 g/L
NaCl	10.0 g/L	5.0 g/L	0 g/L
Total solids	25.0 g/L	20.0 g/L	15.0 g/L
Best for	Standard E. coli — global default	Salt-sensitive strains	Very low osmolarity

Key Features

- Standard Miller formulation — global default for routine E. coli work
- Optimised sodium chloride concentration (10.0 g/L) for osmotic balance
- Supports rapid and reproducible E. coli growth to high optical density
- Suitable as liquid culture medium and as a base for LB Agar preparation
- Easily modified with antibiotics, IPTG, sugars, or other supplements
- Compatible with all standard E. coli K-12 and B strains

Applications

Molecular Biology & Recombinant DNA

- Routine cultivation of E. coli for cloning and plasmid propagation
- Overnight liquid cultures for miniprep, midiprep, and maxiprep
- Preparation of competent cells (chemical or electrocompetent)
- Plasmid maintenance, amplification, and expression
- Recovery broth after transformation (SOC can be substituted)



PRODUCT DESCRIPTION

AS-1270 | LB Broth (Miller)
www.ausamics.com.au

Protein Expression & Fermentation

- Pre-induction liquid culture for recombinant protein expression (BL21, Rosetta)
- Shake flask fermentation for small-scale protein production
- Microbial fermentation and research applications

Teaching & General Laboratory Use

- Standard liquid medium for molecular biology instruction
- General bacterial growth and propagation
- Quality control of competent cell preparations

Preparation Instructions

1. Dissolve 25.0 g of dehydrated medium in 1 litre of demineralised or purified water.
2. Mix or heat gently with agitation until fully dissolved.
3. Verify and adjust pH to 7.0 ± 0.2 if required using 1 M NaOH or 1 M HCl.
4. Sterilise by autoclaving at 121 °C for 15 minutes.
5. Cool to room temperature before use or inoculation.
6. Add filter-sterilised antibiotics or supplements aseptically as required.

Note: To prepare LB Agar (Miller), add 15.0 g/L agar before autoclaving. Cool to 45–50 °C before adding antibiotics and pouring plates.

Common Antibiotic Supplements

Antibiotic	Working Concentration	Stock Concentration	Storage
Ampicillin	100 µg/mL	100 mg/mL in water	-20 °C
Kanamycin	50 µg/mL	50 mg/mL in water	-20 °C
Chloramphenicol	25 µg/mL	25 mg/mL in ethanol	-20 °C
Tetracycline	12.5 µg/mL	12.5 mg/mL in ethanol	-20 °C, protect from light
Gentamicin	10 µg/mL	10 mg/mL in water	-20 °C
Spectinomycin	50 µg/mL	50 mg/mL in water	-20 °C



PRODUCT DESCRIPTION

AS-1270 | LB Broth (Miller)
www.ausamics.com.au

Incubation & Typical Growth

Inoculate with a single colony or from a frozen stock. Incubate with shaking (200–250 rpm) at 37 °C for 12–16 hours for overnight cultures, or 3–4 hours for mid-log phase (OD₆₀₀ 0.4–0.8).

Organism	Expected Result	Incubation
E. coli K-12 (e.g. DH5α, TOP10)	Good growth — turbid broth, OD ₆₀₀ >2.0 overnight	37 °C, shaking, 12–16 h
E. coli BL21(DE3)	Good growth — suitable for protein expression	37 °C, shaking, 3–4 h to OD 0.6
E. coli DH10B	Good growth — high-efficiency transformation host	37 °C, shaking, 12–16 h
Salmonella typhimurium ATCC 14028	Good growth	37 °C, shaking, 12–16 h

Storage & Stability

Dehydrated powder	Store at 15–30 °C in a dry, tightly closed container
Protect from	Moisture, direct light, and extreme temperatures
Prepared broth	2–8 °C, use within 1 week
Shelf life (dehydrated)	As per labelled expiry date
Do not use if	Contamination, precipitation, turbidity, or pH shift observed

Quality & Compliance

Manufactured under controlled conditions for batch-to-batch consistency. Performance tested against ATCC reference organisms prior to release. Compatible with protocols described in Miller (1972) *Experiments in Molecular Genetics*, Sambrook & Russell (*Molecular Cloning*), and *Current Protocols in Molecular Biology*.

Reference

Miller, J.H. (1972). *Experiments in Molecular Genetics*. Cold Spring Harbor Laboratory, New York.



PRODUCT DESCRIPTION

AS-1270 | LB Broth (Miller)
www.ausamics.com.au

Customs & Trade Information

HS / AHECC Code	3821.00.00
Description	Prepared culture media for the development or maintenance of microorganisms
Country of Origin	Australia

Disclaimer

This product is manufactured and supplied by AuSaMicS Pty Ltd for laboratory and research use only. It is not intended for human or veterinary consumption, therapeutic use, or in vitro diagnostic procedures without appropriate validation. AuSaMicS Pty Ltd makes no representations or warranties, express or implied, regarding the fitness of this product for any particular purpose beyond its stated intended use. Users are solely responsible for compliance with all applicable laws, regulations, and safety requirements. AuSaMicS Pty Ltd shall not be liable for any direct, indirect, incidental, or consequential damages arising from the use or misuse of this product. All information provided is believed to be accurate at the time of publication and is subject to change without notice.