



## PRODUCT DESCRIPTION

AS-1338 | Rappaport Vassiliadis Broth (RVS)  
www.ausamics.com.au

### Rappaport-Vassiliadis Soya (RVS) Broth | AS-1338 |

Selective Enrichment for *Salmonella* spp. from Food, Environmental, and Clinical Specimens

**Catalogue Number:** AS-1338 | acc. harm. EP / USP / JP

Target Organism	Pharmacopoeial	Final pH	HS Code
Salmonella spp.	EP / USP / JP harmonised	5.2 ± 0.2 at 25 °C	3821.00.00

#### Overview

**Rappaport Vassiliadis Broth (RVS)** — also designated Rappaport-Vassiliadis Soya Peptone Broth — is a selective liquid enrichment medium formulated for the isolation of *Salmonella* spp. from food, animal feed, environmental samples, and clinical specimens. The medium was originally described by Vassiliadis et al. (1981) as a modification of the earlier Rappaport formula (1954), substituting soy peptone for casein digest to improve recovery of stressed and low-inoculum *Salmonella*.

The medium is harmonised across major international pharmacopoeias — EP 2.6.13, USP <62>, and JP Chapter 4.05 — for Microbial Examination of Non-Sterile Products: Tests for Specified Microorganisms (*Salmonella*). It is also the standard secondary enrichment medium prescribed by ISO 6579-1:2017 (food and animal feed), ISO 19250 (water), and APHA Standard Methods for detection of *Salmonella*.

#### Mode of Action

RVS Broth achieves *Salmonella* selectivity through a combination of four mechanisms: (1) **Malachite green** — a triphenylmethane dye that inhibits Gram-positive organisms and most Gram-negative non-*Salmonella* bacteria at the 0.036 g/L concentration used; (2) **Magnesium chloride** — raises osmotic pressure to ~1400 mOsm/kg, exploiting *Salmonella*'s superior tolerance to elevated  $Mg^{2+}$  and osmotic stress compared to competing flora; (3) **Acidic pH (5.2)** — further inhibits most competing Gram-negative bacteria while *Salmonella* remains viable; (4) **Soy peptone** — provides a balanced amino acid and growth factor profile that preferentially supports *Salmonella* enrichment.

#### Typical Composition (per litre)

Ingredient	CAS Number	Function	Amount (g/L)
Soy Peptone (Tryptone Soya Peptone)	73049-73-7	Sole nitrogen and carbon source — balanced amino acids for <i>Salmonella</i>	5.0



## PRODUCT DESCRIPTION

AS-1338 | Rappaport Vassiliadis Broth (RVS)  
www.ausamics.com.au

Sodium Chloride (NaCl)	7647-14-5	Base osmotic balance	8.0
Potassium Dihydrogen Phosphate (KH <sub>2</sub> PO <sub>4</sub> )	7778-77-0	pH buffer (acidic component)	1.6
Dipotassium Hydrogen Phosphate (K <sub>2</sub> HPO <sub>4</sub> )	7758-11-4	pH buffer (alkaline component)	0.8
Magnesium Chloride Hexahydrate (MgCl <sub>2</sub> ·6H <sub>2</sub> O)	7791-18-6	Elevated osmotic pressure — selective for Salmonella	13.4
Malachite Green Oxalate	2437-29-8	Selective inhibitor — Gram-positive and non-Salmonella Gram-negatives	0.036
Purified Water	7732-18-5	Solvent	q.s. 1 L

**Total (without water):** 28.836 g/L | Final pH: 5.2 ± 0.2 at 25 °C

### Key Features

- Pharmacopoeial harmonised — acc. EP 2.6.13, USP <62>, JP Chapter 4.05
- ISO 6579-1:2017 compliant — international standard for Salmonella in food and animal feed
- Dual selectivity mechanism — malachite green + elevated osmotic pressure (MgCl<sub>2</sub>) + acidic pH
- Soy peptone base — superior Salmonella recovery compared to casein peptone variants
- Supports recovery of stressed, freeze-damaged, and heat-injured Salmonella cells
- Used as secondary selective enrichment following non-selective pre-enrichment (BPW)
- Applicable to food, animal feed, water, environmental swabs, and pharmaceutical non-sterile products

### Applications

#### Food & Animal Feed Safety

- Secondary selective enrichment of Salmonella from pre-enriched BPW cultures — ISO 6579-1:2017
- Broad food matrices: raw meat, poultry, eggs, dairy, dried foods, spices, herbs, pet food
- Animal feed and raw material Salmonella screening per ISO 6579-1
- Environmental swab enrichment from food production facilities

#### Pharmaceutical Quality Control

- Salmonella testing of non-sterile pharmaceutical products — EP 2.6.13, USP <62>, JP Chapter 4.05
- Testing of herbal medicines, excipients, and raw botanical materials
- Finished product release testing for Salmonella absence per pharmacopoeial harmonised method



## PRODUCT DESCRIPTION

AS-1338 | Rappaport Vassiliadis Broth (RVS)  
www.ausamics.com.au

### Water & Environmental Testing

- Salmonella enrichment from drinking water, wastewater, surface water — ISO 19250
- Environmental monitoring swabs from processing and manufacturing facilities
- Sewage sludge and biosolids Salmonella screening

### Workflow — Salmonella Detection (ISO 6579-1)

Step	Medium	Incubation	Purpose
1. Non-selective pre-enrichment	Buffered Peptone Water (BPW)	37 °C, 18 ± 2 h	Resuscitate stressed/injured Salmonella cells
2. Selective enrichment	RVS Broth (AS-1338) ← this product	41.5 °C, 24 ± 3 h	Select and amplify Salmonella; suppress competing flora
3. Selective plating	XLD Agar + Chromogenic Salmonella Agar	37 °C, 24–48 h	Presumptive colony identification
4. Confirmation	Biochemical + serological	—	Confirmed positive result

*Note: Transfer 0.1 mL of BPW pre-enrichment to 10 mL RVS Broth (1:100 inoculation ratio as specified in ISO 6579-1:2017).*

### Preparation Instructions

1. Dissolve 26.75 g (or as stated on label) of dehydrated RVS Broth in 1 litre of purified water.
2. Mix with gentle agitation at room temperature — do NOT heat. Malachite green and MgCl<sub>2</sub> are heat-sensitive.
3. Dispense into appropriate volumes: 10 mL per tube (for 0.1 mL inoculum) or 100 mL per flask.
4. Sterilise by autoclaving at 115 °C for 15 minutes. Do NOT autoclave at 121 °C — excessive heat degrades malachite green and alters osmolarity.
5. Cool to room temperature. Medium should appear bluish-green and clear.
6. Store at 2–8 °C, protected from light. Use within 3 months.

*Critical: Autoclave at 115 °C (not 121 °C). Higher temperature irreversibly degrades malachite green, reducing selectivity and potentially causing false-negative results.*

### Storage & Stability



## PRODUCT DESCRIPTION

AS-1338 | Rappaport Vassiliadis Broth (RVS)  
www.ausamics.com.au

<b>Dehydrated powder</b>	15–30 °C, tightly sealed, dry, protected from light
<b>Prepared medium</b>	2–8 °C, protected from light; use within 3 months
<b>Shelf life (powder)</b>	As per labelled expiry date
<b>Do not use if</b>	Colour changed (should be bluish-green), turbid, or precipitate visible before inoculation

## Customs & Trade Information

<b>HS / AHECC Code</b>	3821.00.00
<b>Description</b>	Prepared culture media for the development or maintenance of microorganisms
<b>Country of Origin</b>	Australia

## Literature & References

- Vassiliadis, P., Kalapothaki, V., Trichopoulos, D., Papadakis, J.A. & Serie, C. (1981). Salmonella isolation with Rappaport's Vassiliadis medium. J. Appl. Bacteriol., 51(2), 233–239.
- Rappaport, F., Konforti, N. & Navon, B. (1954). A new enrichment medium for certain Salmonellae. J. Clin. Pathol., 9(3), 261–266.
- ISO 6579-1:2017. Microbiology of the food chain — Horizontal method for the detection, enumeration and serotyping of Salmonella — Part 1: Detection. ISO, Geneva.
- ISO 19250:2010. Water quality — Detection of Salmonella spp. ISO, Geneva.
- European Pharmacopoeia, Chapter 2.6.13 — Microbial examination of non-sterile products: Tests for specified microorganisms.
- USP <62> — Microbiological examination of non-sterile products: Tests for specified microorganisms.
- JP Chapter 4.05 — Microbial Limit Test II.
- APHA Standard Methods for the Examination of Water and Wastewater (2017), Method 9260 B.

## Disclaimer

*This product is manufactured and supplied by AuSaMics Pty Ltd for laboratory and research use only. It is not intended for human or veterinary consumption, therapeutic use, or in vitro diagnostic procedures without appropriate validation. AuSaMics Pty Ltd makes no representations or warranties, express or implied, regarding the fitness of this product for any particular purpose beyond its stated intended use. Users are solely responsible for compliance with all applicable laws, regulations, and safety requirements. AuSaMics Pty Ltd shall not be liable for any direct, indirect, incidental, or consequential damages arising from the use or misuse of this product. All information provided is believed to be accurate at the time of publication and is subject to change without notice.*