



Technical Data Sheet

TDS-AS-1249
www.ausamics.com.au

Product Identification

Product Name	Helicobacter pylori Selective Medium Base
Catalog No.	AS-1249
Lot No.	HPS260301
Mfg. Date	March 2026
Retest Date	March 2028
Grade	Microbiological / Selective Grade
Base Medium	Columbia Blood Agar Base + selective antibiotic supplement + defibrinated sheep blood

Physical & Chemical Properties

Appearance (powder)	Homogeneous, free-flowing powder; light beige to pale pink
Appearance (prepared)	Opaque, red-brown gel with 5% sheep blood; homogeneous
Odour	Faint, characteristic (blood agar)
pH (prepared, 25 °C)	7.3 ± 0.2
Agar gel / melt	Gels ~32–34 °C / Melts ~84–86 °C
Loss on Drying	≤ 7.0% w/w (105 °C / 2 h)
Haemolysis Pattern	Gamma (non-haemolytic) for H. pylori colonies

Base Medium Composition per Litre (Columbia Base)

Component	Function	Amount (g/L)
Peptone (mixed — casein/meat/heart)	Nitrogen, amino acids — supports fastidious organisms	23.00
Cornstarch (soluble)	Carbon source + protective agent for organism viability	1.00



Technical Data Sheet

TDS-AS-1249
www.ausamics.com.au

Component	Function	Amount (g/L)
Sodium Chloride	Osmotic balance	5.00
Agar (bacteriological grade)	Solidifying agent	10.00
TOTAL (base powder per litre)		39.00
Defibrinated sheep blood	Growth factors (haemin, NAD), haemolysis indicator — SUPPLEMENT	5–10% v/v

Antibiotic Selective Supplement

Antibiotic	Concentration (typical)	Target Organisms
Vancomycin	10 mg/L	Gram-positive bacteria
Trimethoprim	5 mg/L	Enteric Gram-negative bacilli
Cefsulodin	5 mg/L	<i>Pseudomonas aeruginosa</i>
Amphotericin B	2 mg/L	Yeasts and moulds

Preparation Protocol

Step 1 — Suspension	Suspend 39 g of dehydrated Columbia base in 1 L of purified/distilled water.
Step 2 — Dissolution	Heat with agitation until completely dissolved. Bring to a rolling boil for 1 minute.
Step 3 — Sterilisation	Autoclave at 121 °C (15 psi) for 15 minutes. Do not over-autoclave.
Step 4 — Cooling	Cool to 45–50 °C in a water bath. Temperature is critical — above 50 °C will lyse blood cells.
Step 5 — Blood addition	Aseptically add 5–10% v/v defibrinated sheep blood (pre-warmed to 45 °C). Mix gently — avoid frothing.
Step 6 — Antibiotics	Add antibiotic selective supplement aseptically per manufacturer's instructions. Mix gently.
Step 7 — Pouring	Pour 15–20 mL per 90 mm Petri dish under laminar flow. Allow to solidify on a level surface.



Technical Data Sheet

TDS-AS-1249
www.ausamics.com.au

Step 8 — Drying

Dry plates at 37 °C for 30–45 min. Store at 2–8 °C in sealed bags. Use within 2 weeks.

Incubation Conditions

Temperature	37 °C (optimal); up to 42 °C for primary clinical isolates
Atmosphere	Microaerophilic: 5–10% CO ₂ , 5–10% O ₂ , balance N ₂ (CampyPak or equivalent generator)
Duration	3–7 days; examine at 72 h then every 24 h
Humidity	≥ 95% RH — critical; prevent plate desiccation by sealing or humidified incubation
Do NOT use	Aerobic or anaerobic conditions — <i>H. pylori</i> will not grow

Quality Control Specifications

Parameter	Specification	Method
Appearance (powder)	Homogeneous; light beige to pale pink; no lumps	Visual inspection
pH (prepared, 25 °C)	7.1 – 7.5	Potentiometry (ISO 10523)
Loss on Drying	≤ 7.0% w/w	105 °C / 2 h (gravimetric)
Growth — <i>H. pylori</i> ATCC 43504	Good growth; small translucent colonies ≤ 5 days	Microaerophilic culture
Inhibition — <i>E. coli</i> ATCC 25922	Complete / ≥ 99% inhibition	Selectivity test
Inhibition — <i>Candida albicans</i> ATCC 10231	Complete inhibition	Antifungal selectivity
Haemolysis pattern	Gamma (non-haemolytic) for <i>H. pylori</i> colonies	Visual — plate check
Sterility (prepared medium)	No growth at 14 days (30–35 °C, aerobic)	Incubation sterility check

Technical References



Technical Data Sheet

TDS-AS-1249
www.ausamics.com.au

Primary Reference	Marshall & Warren (1984). Lancet, 323:1311–1315
Clinical Standard	Maastricht V/Florence Consensus (2017). Gut, 66(1):6–30
Diagnostic Method	Culture — gold standard for H. pylori AST and genotyping
Safety Reference	IARC Monograph Vol. 61 (1994): H. pylori classified Group I carcinogen
Containment	PC2 laboratory required. Handle per AS/NZS 2243.3:2010 Biosafety in Laboratories

Disclaimer: AuSaMicS Pty Ltd (ABN 56 676 640 467) warrants that this product meets the specifications stated herein at time of manufacture and release. This product is for laboratory and research use only. Not intended for human, animal, food, feed, or pharmaceutical use. AuSaMicS Pty Ltd shall not be held liable for any direct, indirect, incidental, or consequential damages resulting from use outside its intended application. By using this product the purchaser agrees to these terms.