



Technical Data Sheet

TDS-AS-1337
www.ausamics.com.au

Product Identification

| | |
|--------------|---|
| Product Name | R2A Agar (Reasoner's 2A) |
| Catalog No. | AS-1337 |
| Lot No. | R2A260301 |
| Mfg. Date | March 2026 |
| Retest Date | March 2028 |
| Grade | Microbiological / Analytical Grade |
| Standard | ISO 6222:1999 · Ph. Eur. 11th Ed. · APHA SM-9215B |

Physical & Chemical Properties

| | |
|------------------------|--|
| Appearance (powder) | Homogeneous, free-flowing powder; pale cream to light beige |
| Appearance (prepared) | Clear to slightly opalescent gel; light amber colour |
| Odour | Faint, characteristic (peptone/yeast extract) |
| pH (prepared, 25 °C) | 7.2 ± 0.2 |
| Agar gel / melt point | Gels ≈ 32–34 °C / Melts ≈ 84–86 °C |
| Loss on Drying (spec.) | ≤ 7.0% w/w (105 °C / 2 h) |
| Solubility | 18.12 g/L in purified water on heating; swells in cold water |

Composition per Litre — ISO 6222 / Ph. Eur. Standard

| Component | Function | g / L |
|---|--|-------|
| Yeast Extract | Vitamins, amino acids — primary C/N source | 0.50 |
| Proteose Peptone No. 3 (or Polypeptone) | Peptides — slow-release nitrogen source | 0.50 |

| Component | Function | g / L |
|---------------------------------------|--|-------|
| Casein Hydrolysate (acid) | Amino acids — supplementary nitrogen | 0.50 |
| Glucose (Dextrose) | Fermentable carbon source | 0.50 |
| Starch (Soluble) | Carbon source + peroxide neutralisation — protects injured cells | 0.50 |
| K ₂ HPO ₄ | pH buffer — stabilises at 7.2 | 0.30 |
| Sodium Pyruvate | Peroxide scavenger — rescues chlorine-injured cells | 0.30 |
| MgSO ₄ · 7H ₂ O | Divalent cation — enzyme cofactor, membrane integrity | 0.024 |
| Agar (bacteriological grade) | Solidifying agent | 15.00 |
| TOTAL (per litre) | | 18.12 |

Preparation Protocol

| | |
|------------------------|--|
| Step 1 — Suspension | Weigh 18.12 g of dehydrated R2A Agar. Suspend in 1 L of purified/distilled water (type 1 or equivalent). |
| Step 2 — Dissolution | Heat with frequent agitation. Bring to a rolling boil for 1 minute until completely dissolved. Do not over-heat. |
| Step 3 — Sterilisation | Autoclave at 121 °C (15 psi / 103 kPa) for exactly 15 minutes. Over-autoclaving degrades pyruvate and starch. |
| Step 4 — Pouring | Cool in a 45–50 °C water bath. Pour 15–20 mL per 90 mm Petri dish under sterile conditions. |
| Step 5 — Drying | Allow to solidify on a level surface. Dry at 37 °C for 30–45 min before use. Store at 2–8 °C. |

Incubation & Counting Specifications



Technical Data Sheet

TDS-AS-1337
www.ausamics.com.au

| Protocol | Temperature | Duration | Standard |
|----------------------|-------------|-------------|--------------------------------|
| Standard (preferred) | 20–28 °C | 5–7 days | ISO 6222 / Ph. Eur. 2.6.12 |
| Accelerated | 35–37 °C | 44–72 hours | When shorter protocol required |

Count all visible colonies regardless of size, pigmentation, or morphology. Report results as CFU/100 mL or CFU/mL.

Quality Control Specifications

| Parameter | Specification | Method |
|---|---------------------------------------|----------------------------|
| Appearance (powder) | Homogeneous; cream to beige; no lumps | Visual inspection |
| pH (prepared, 25 °C) | 7.0 – 7.4 | Potentiometry (ISO 10523) |
| Loss on Drying | ≤ 7.0% w/w | 105 °C / 2 h (gravimetric) |
| Growth — <i>P. aeruginosa</i> ATCC 9027 | Recovery ≥ 70% vs. reference medium | ISO 6222 growth promotion |
| Growth — <i>S. aureus</i> ATCC 6538 | Recovery ≥ 70% vs. reference medium | ISO 6222 growth promotion |
| Growth — <i>M. luteus</i> ATCC 9341 | Visible colonies ≤ 48 h at 28 °C | Ph. Eur. 2.6.12 |
| Sterility (prepared medium) | No growth at 14 days (30–35 °C) | Incubation sterility check |

Technical References

| | |
|-------------------|--|
| Primary Reference | Reasoner D.J. & Geldreich E.E. (1985). Appl. Environ. Microbiol. 49(1):1–7 |
| ISO Standard | ISO 6222:1999 — Enumeration of culturable micro-organisms |
| Pharmacopoeial | Ph. Eur. 11th Ed. 2.6.12 — Medium S |
| APHA/AWWA | Standard Methods for Examination of Water & Wastewater, 9215 B |
| US EPA | EPA Method 1604 (selected applications) |



Technical Data Sheet

TDS-AS-1337

www.ausamics.com.au

Disclaimer: *AuSaMicS Pty Ltd (ABN 56 676 640 467) warrants that this product meets the specifications stated herein at time of manufacture and release. This product is for laboratory and research use only. AuSaMicS Pty Ltd shall not be held liable for any direct, indirect, incidental, or consequential damages resulting from use outside its intended application. By using this product the purchaser agrees to these terms.*